

CompactDry Interpretation Tips




Copyright © Shimadzu Diagnostics Corporation All rights reserved.


CompactDry Interpretation Tips

#	Colony Characteristics
1	Food Residual
2	Spread Colony
3	Whole Surface Colored
4	Mass Colonies
5	Mixed Color on EC
6	Light Colonies
7	SL Interpretations
8	White Colonies
9	Different Colors' Colonies


Food Residual

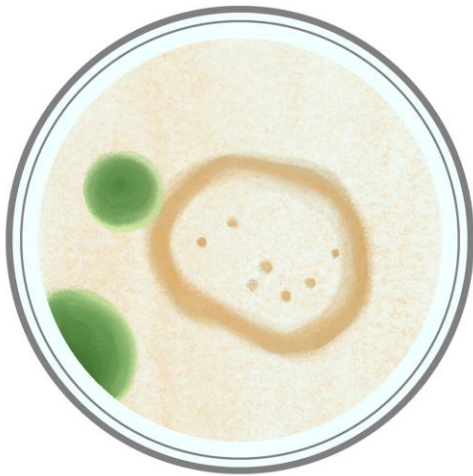

EC

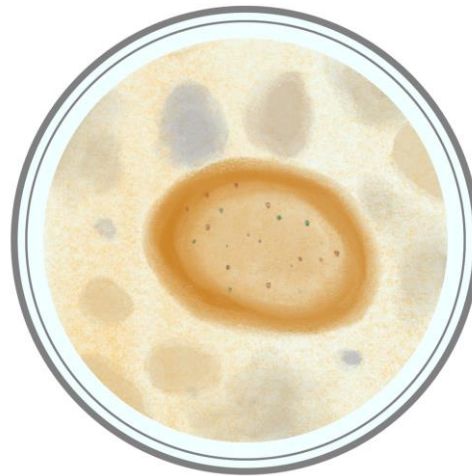


XSA

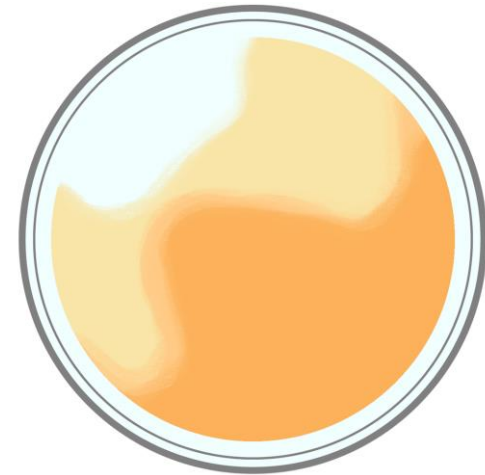


YMR


Food Residual

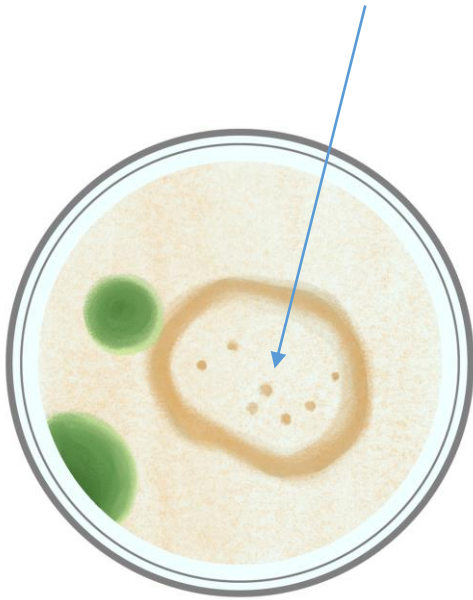

EC

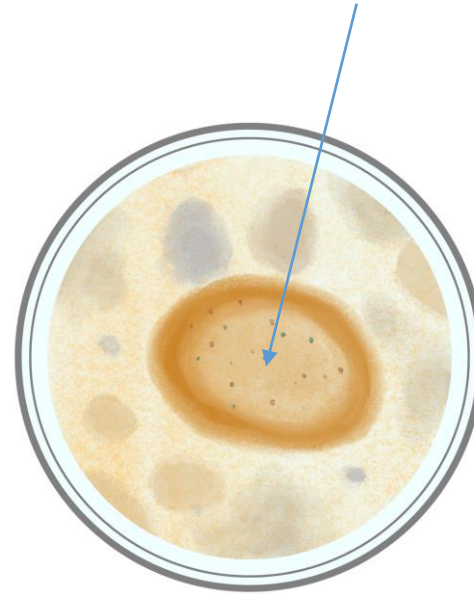


XSA

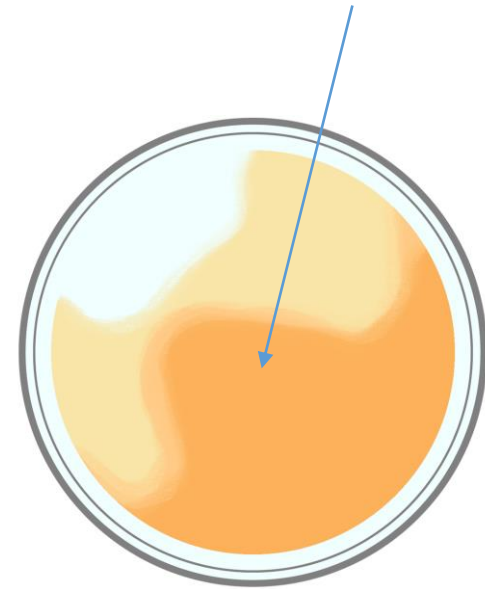


YMR

Spread Colony

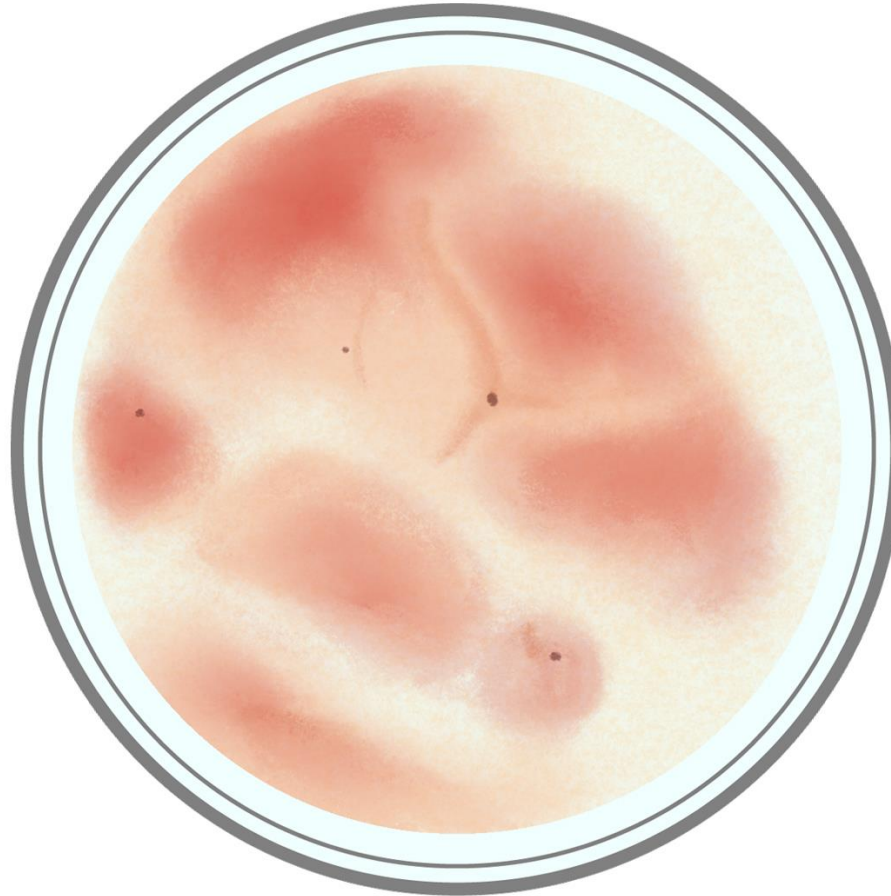


TC

Spread Colony

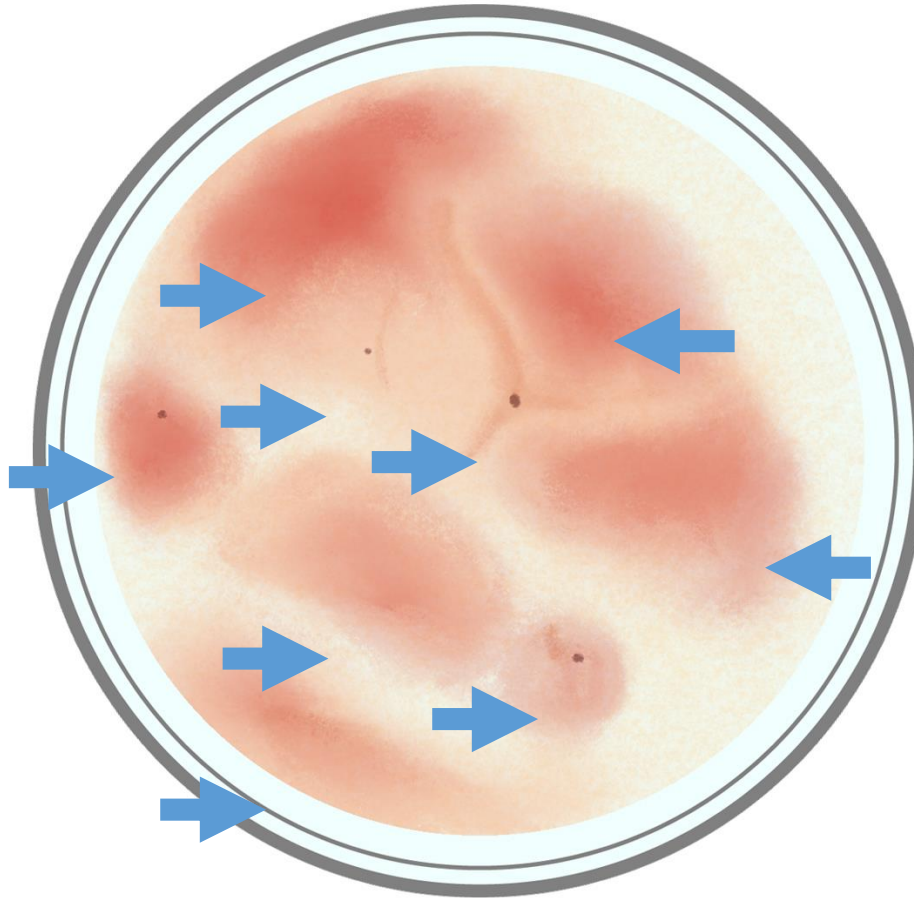


TC

Spread Colony

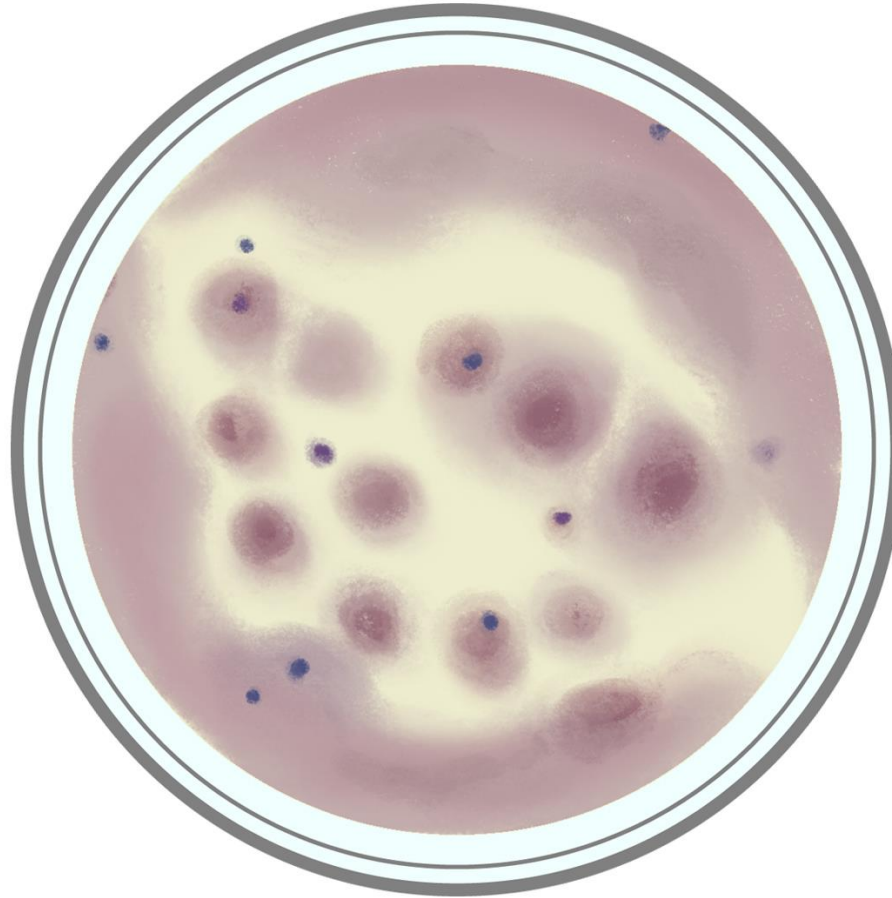


EC

Spread Colony

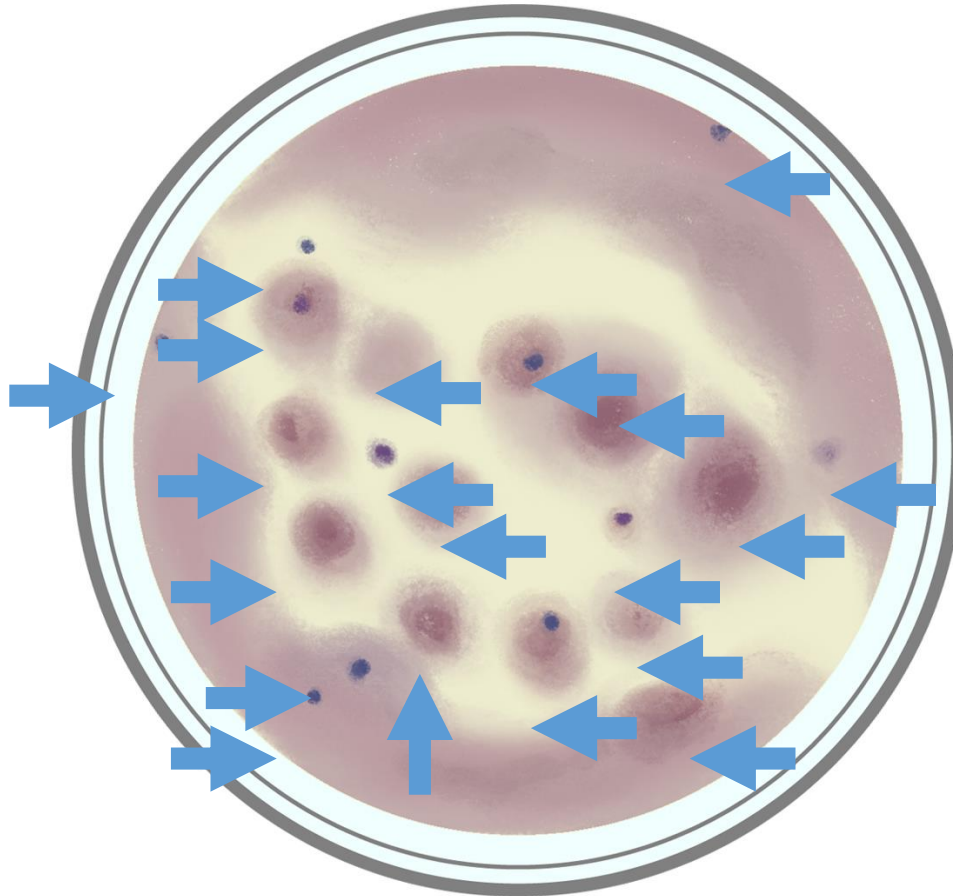


EC

Spread Colony

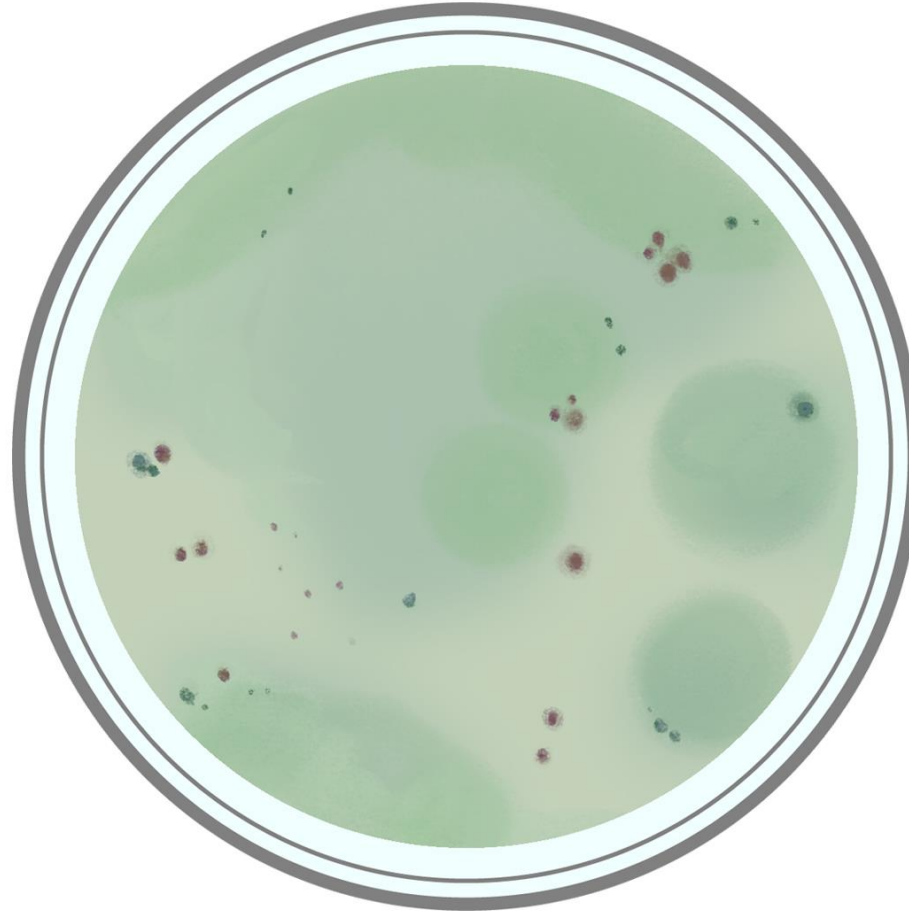


VP

Spread Colony

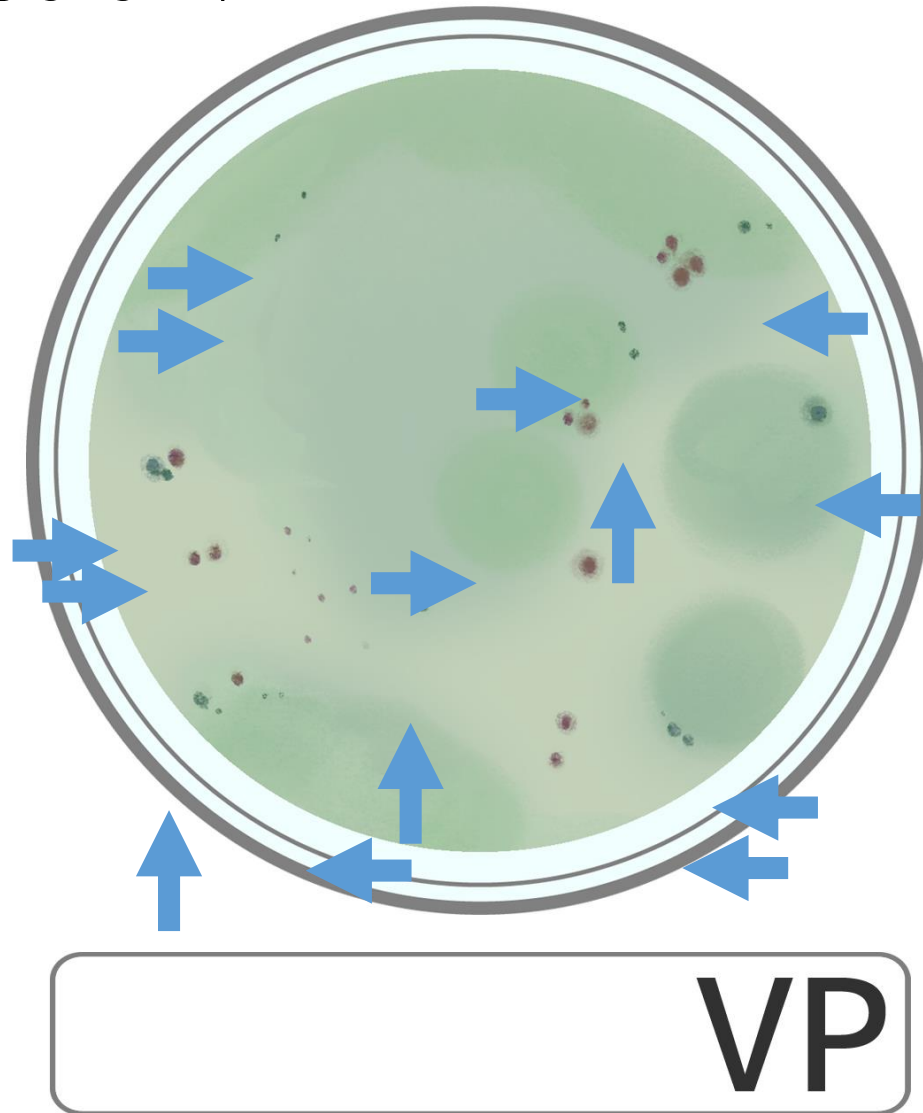


Whole Surface Colored

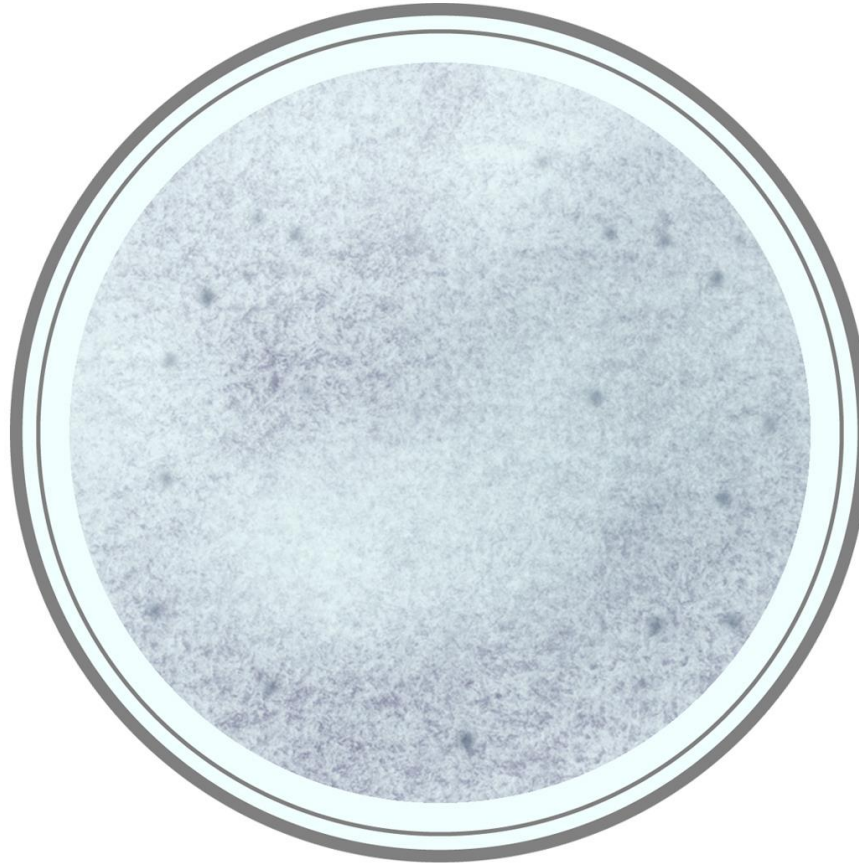


EC

Whole Surface Colored

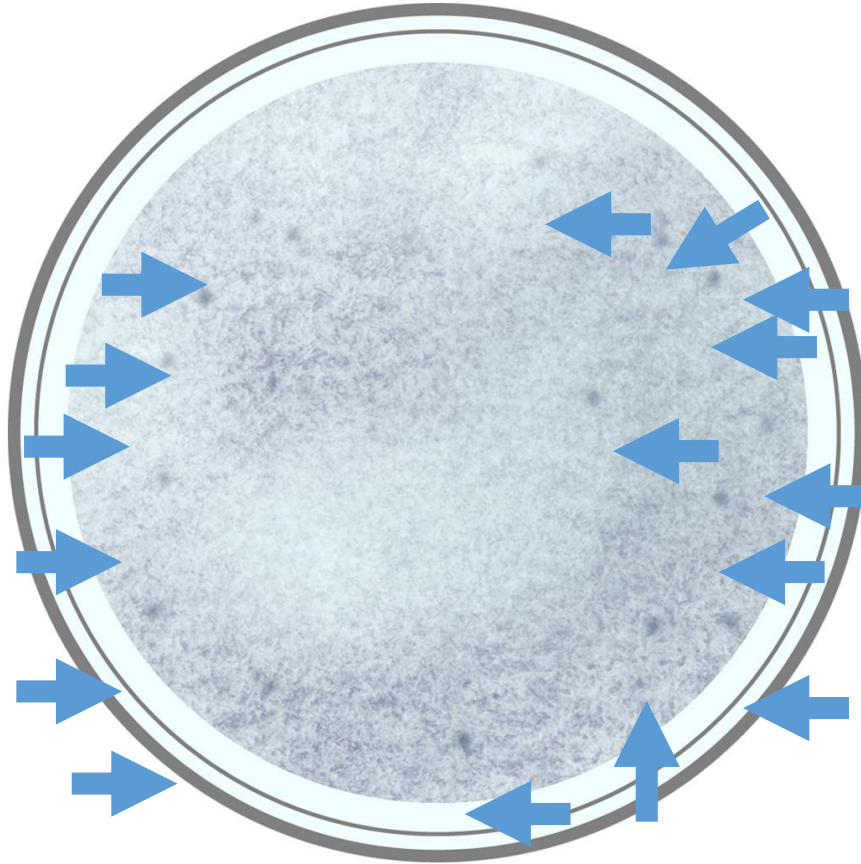


EC

Whole Surface Colored

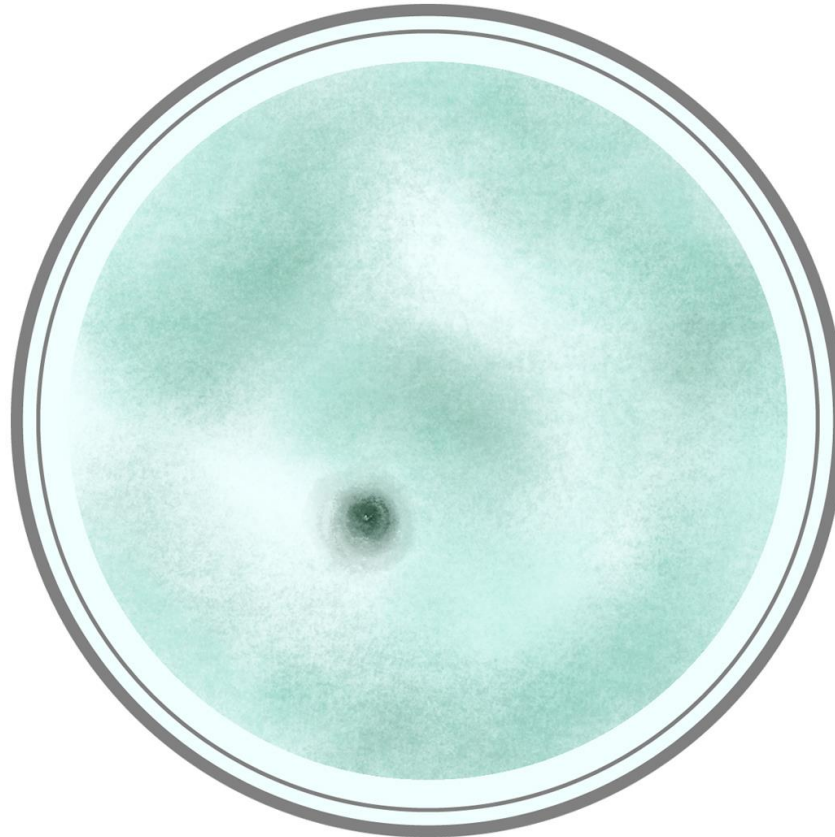


YM

Whole Surface Colored

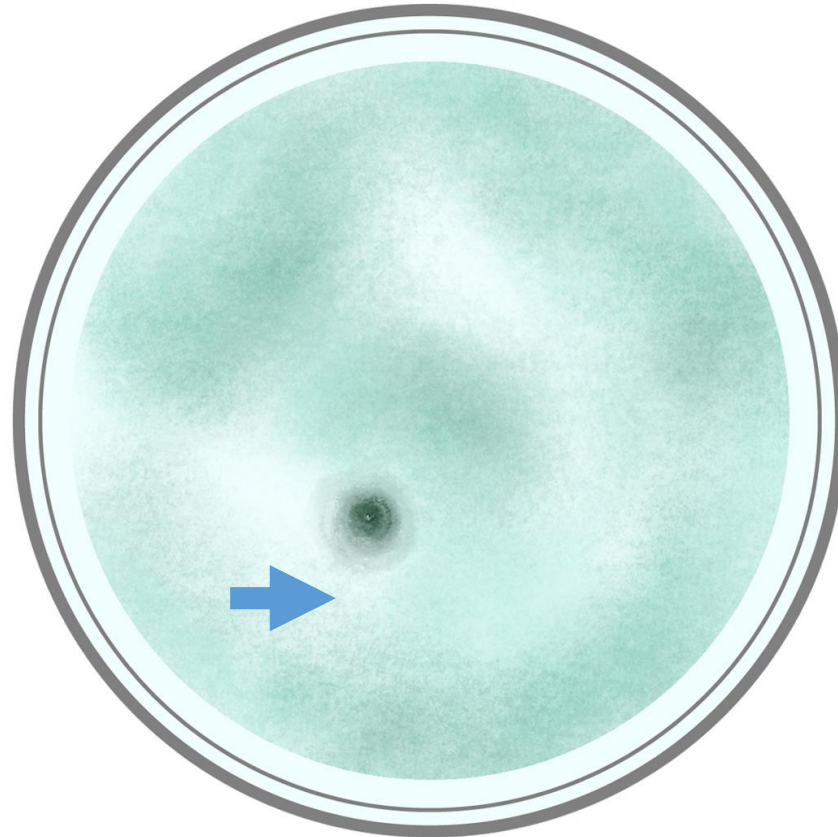


YM

Whole Surface Colored

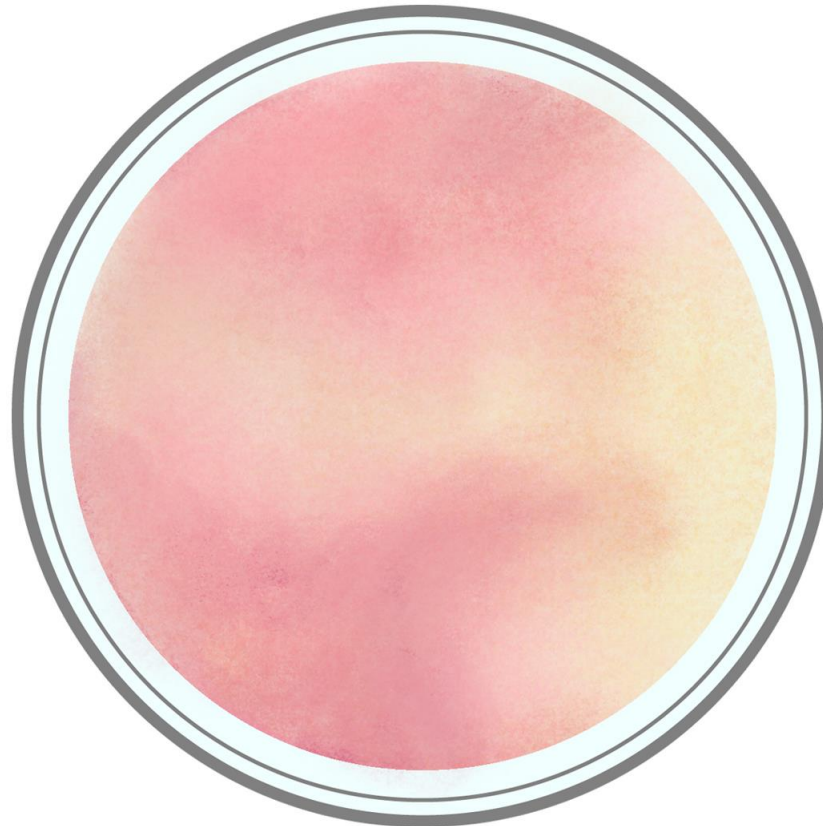


ETB

Whole Surface Colored

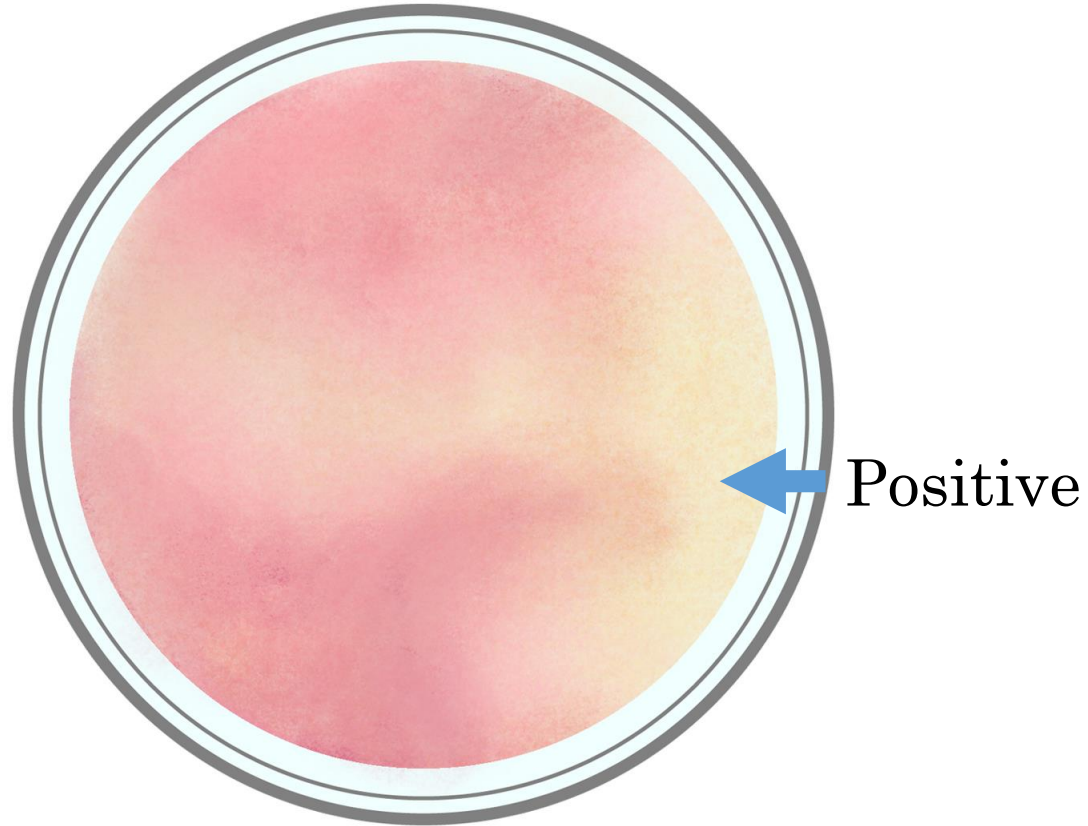


ETB

Mass Colonies

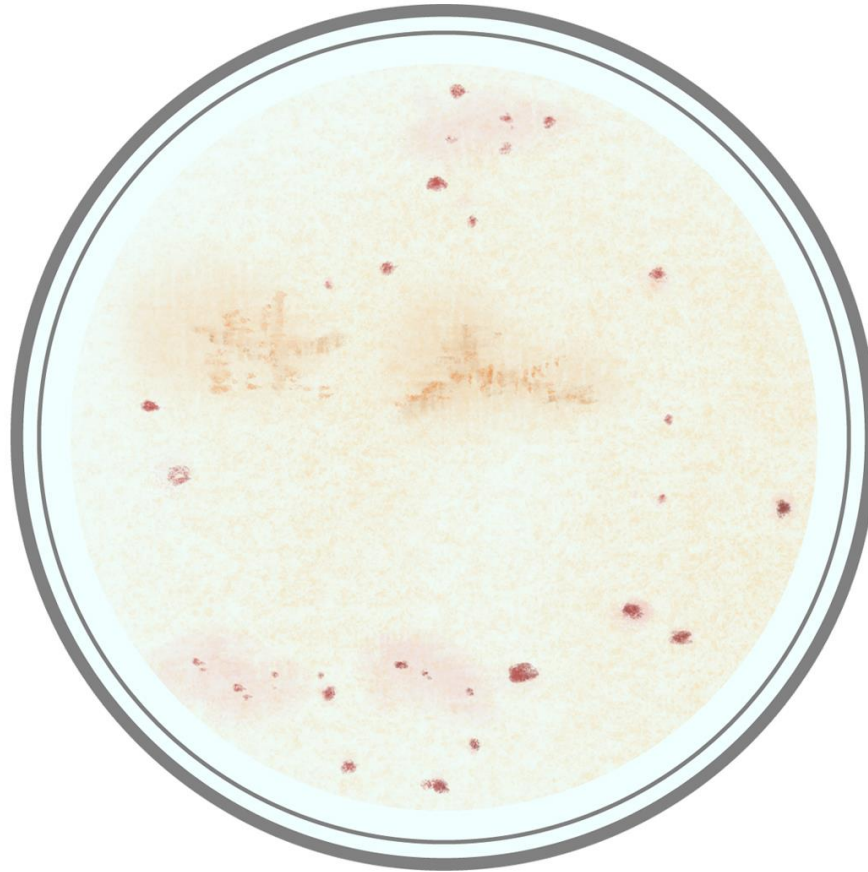


TC

Mass Colonies

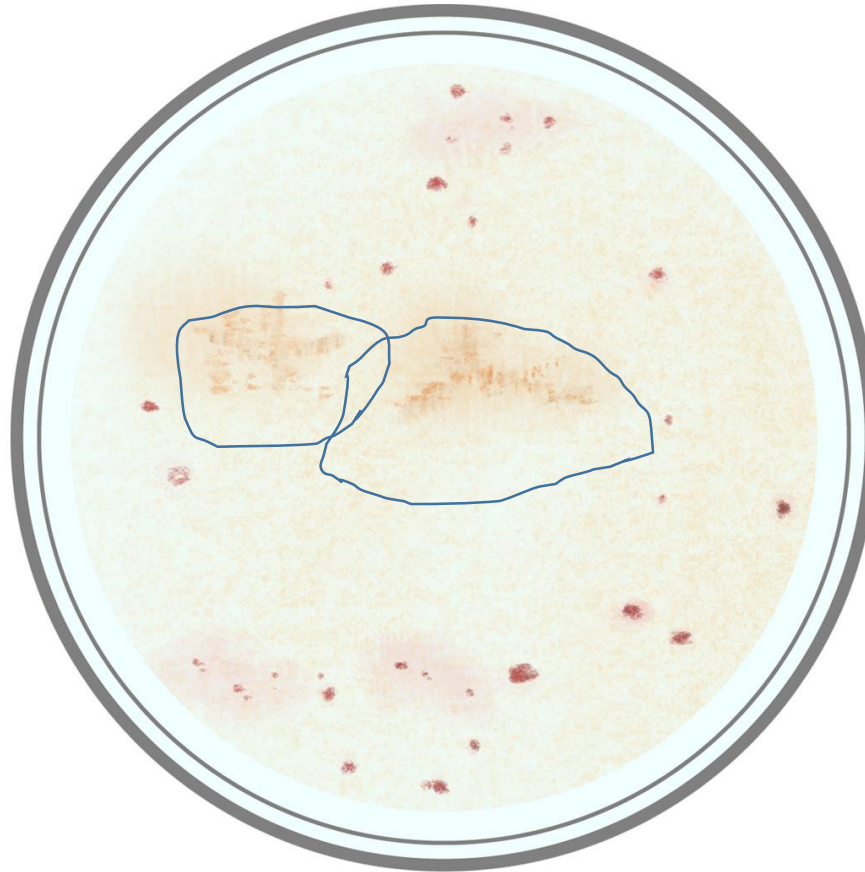


TC

Mass Colonies

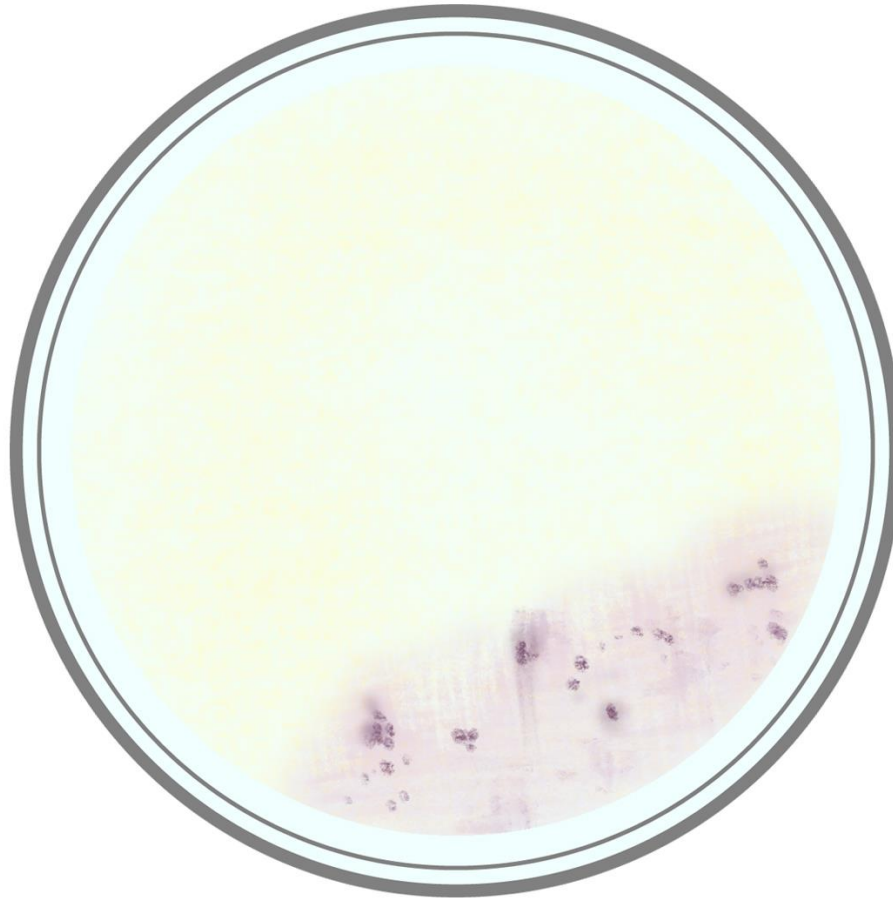


EC

Mass Colonies

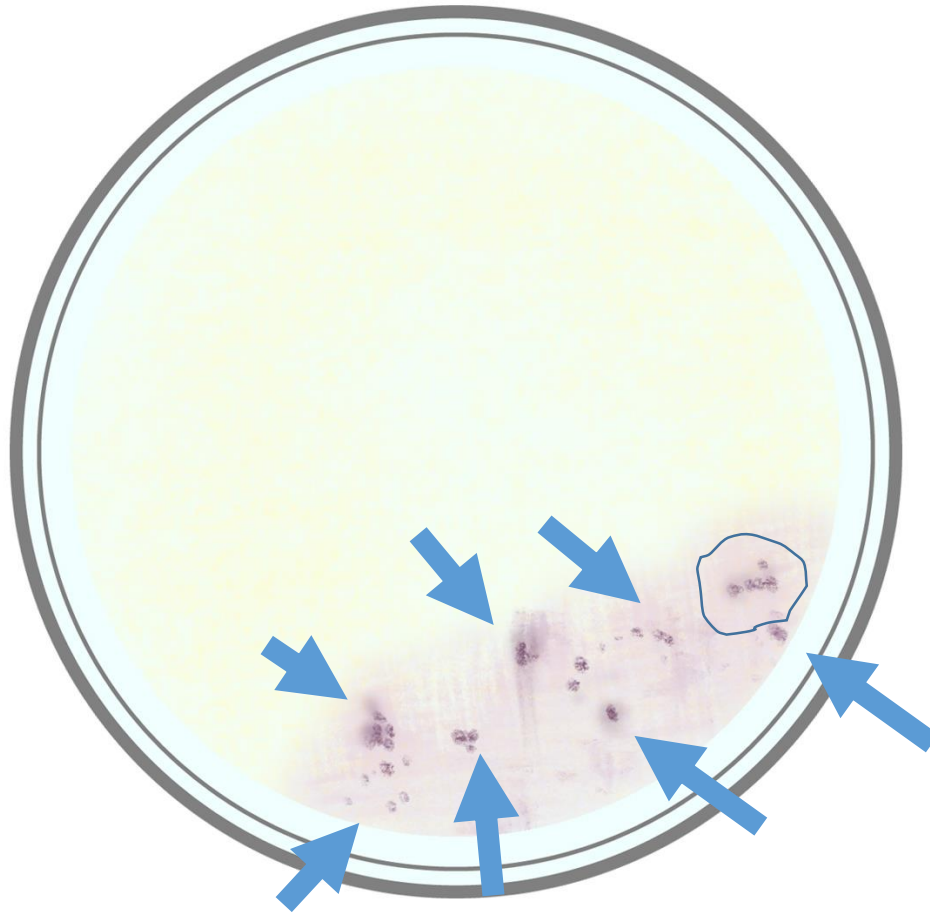


EC

Mass Colonies

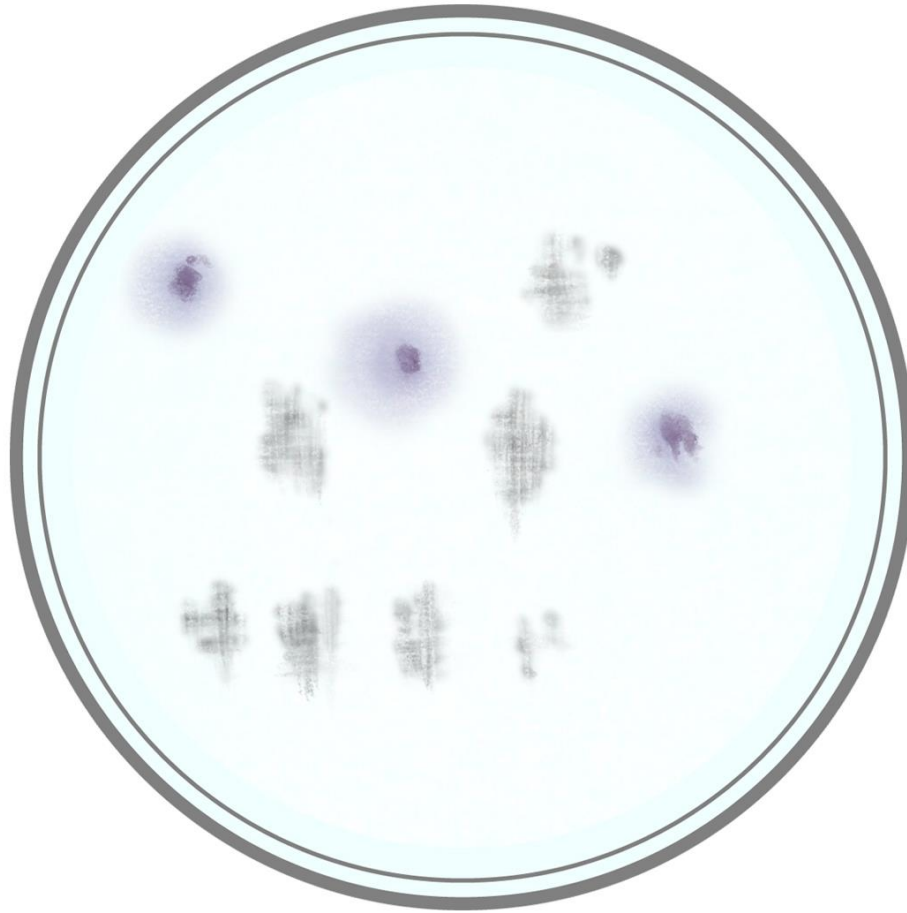


EC

Mass Colonies

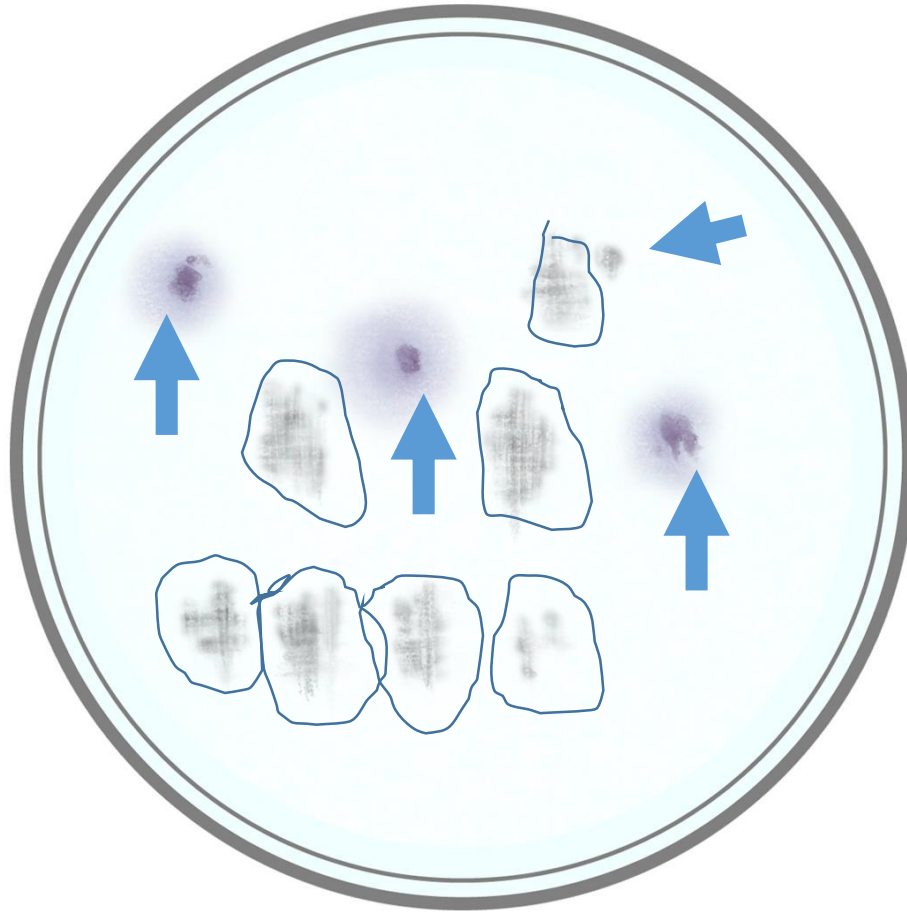


EC


Mixed Color on EC

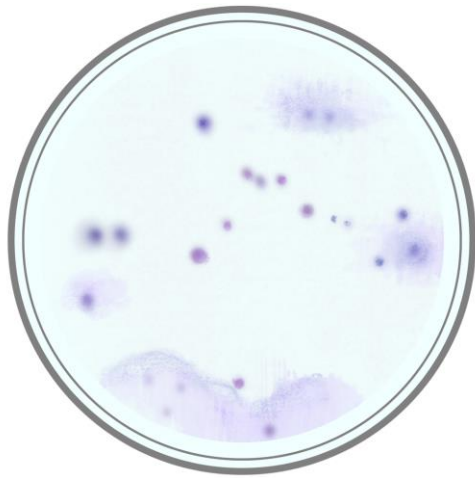

EC

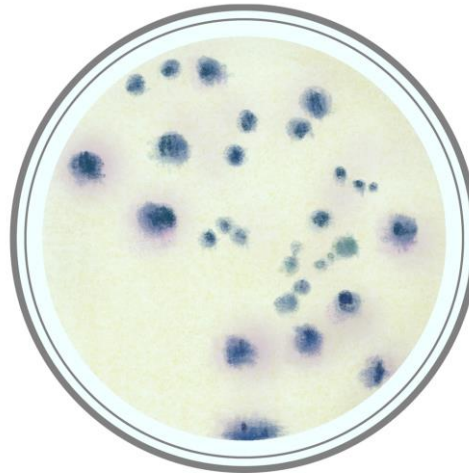


EC

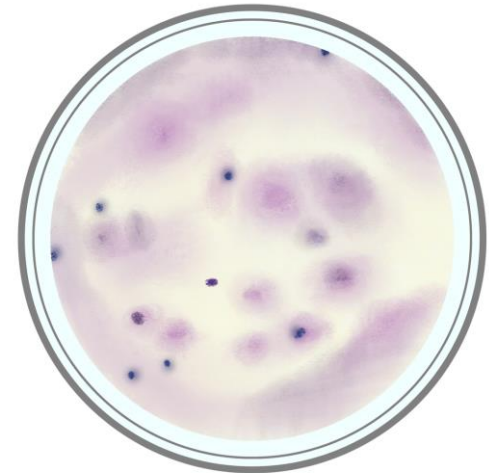


EC

Mixed Color on EC

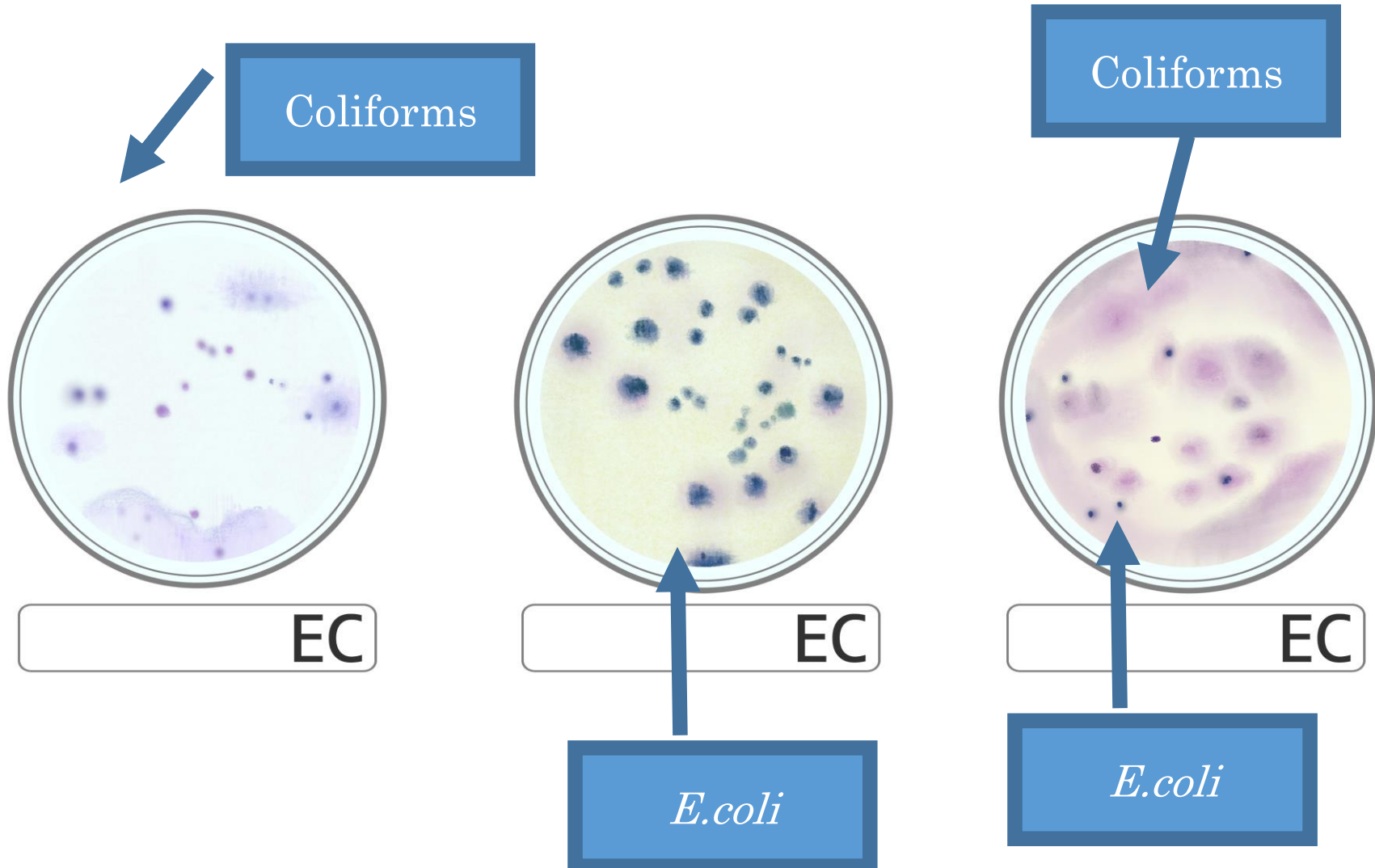


Light Colonies

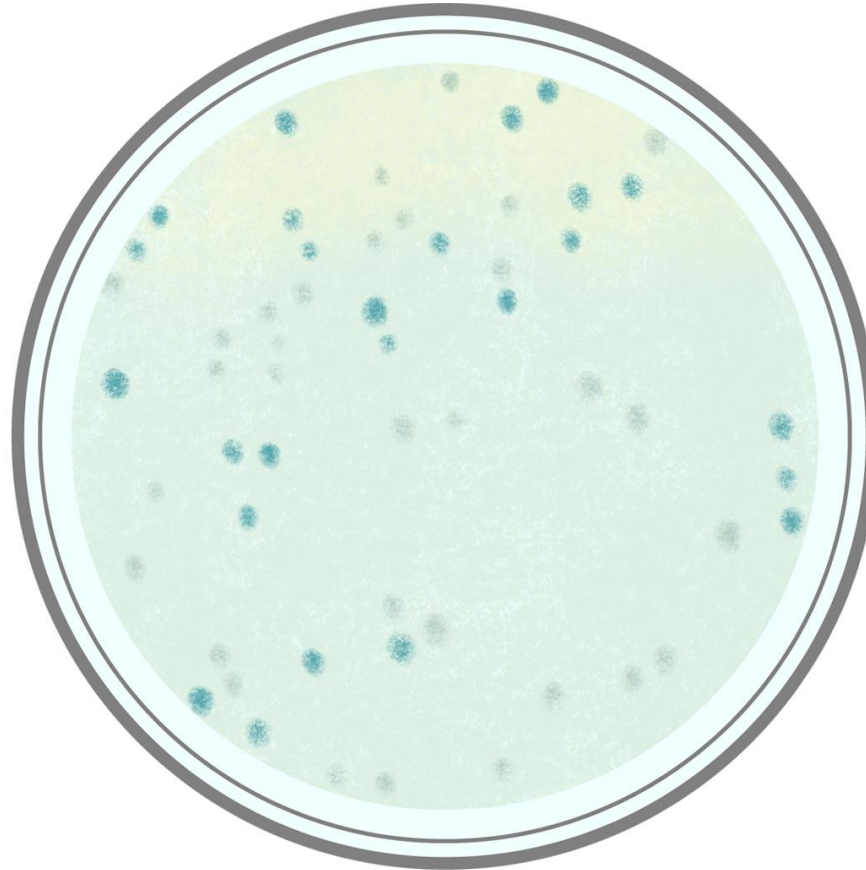

XSA

Light Colonies

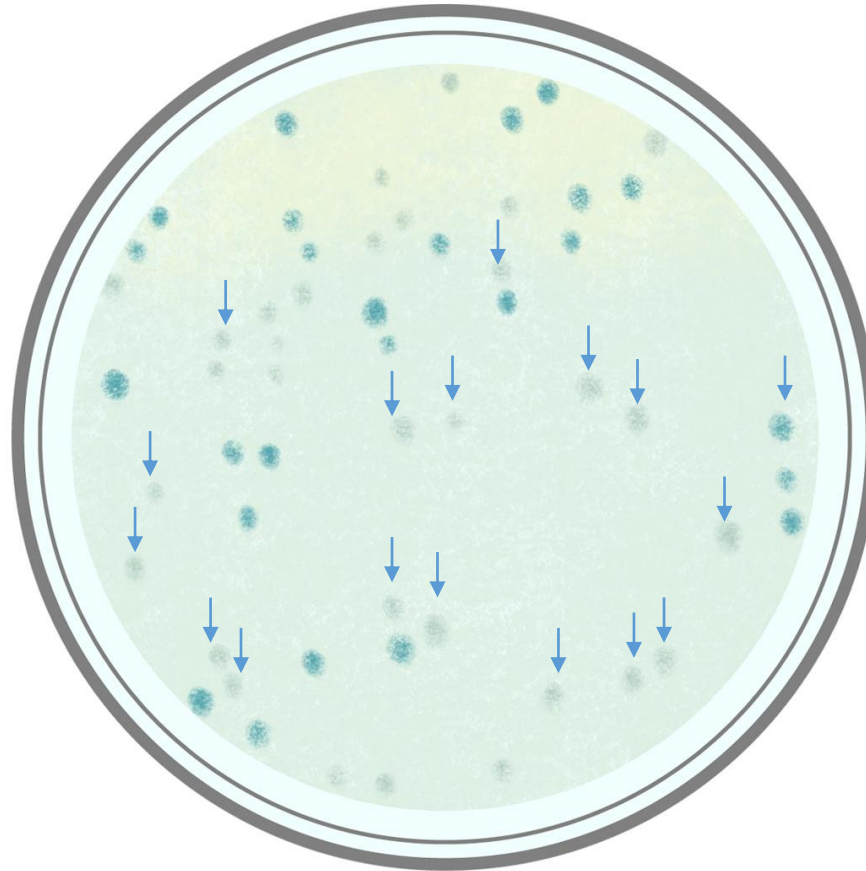


XSA

Light Colonies




LS

Light Colonies

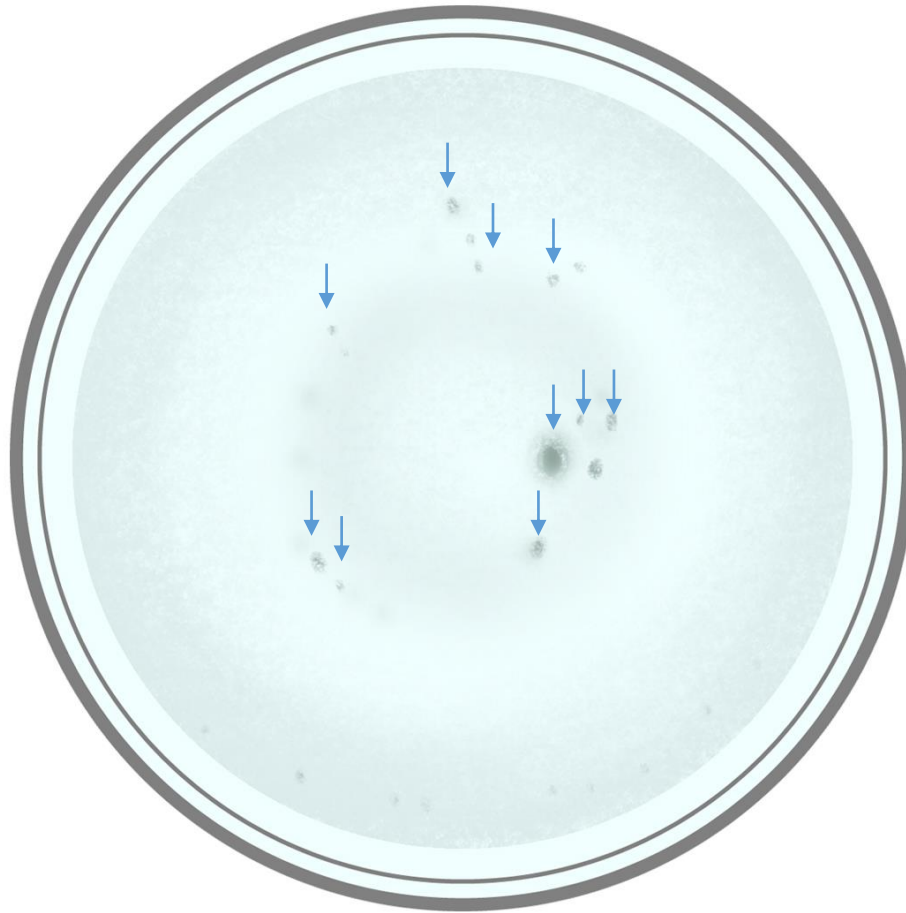


LS

Light Colonies

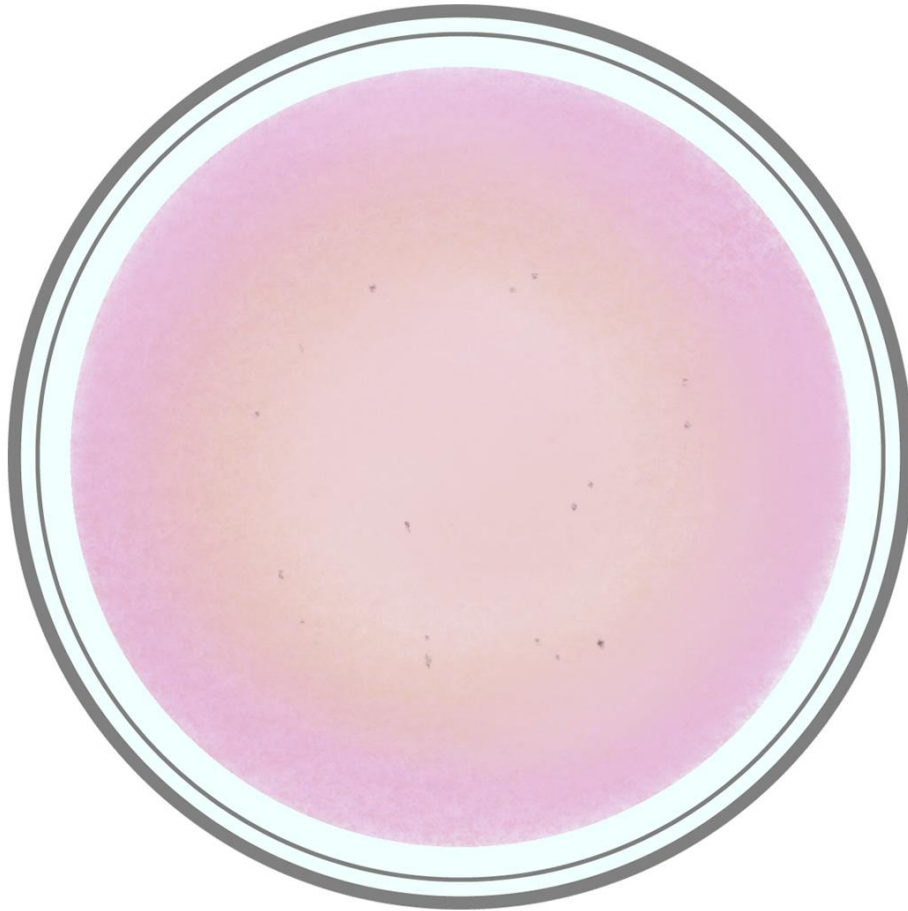

ETB

Light Colonies

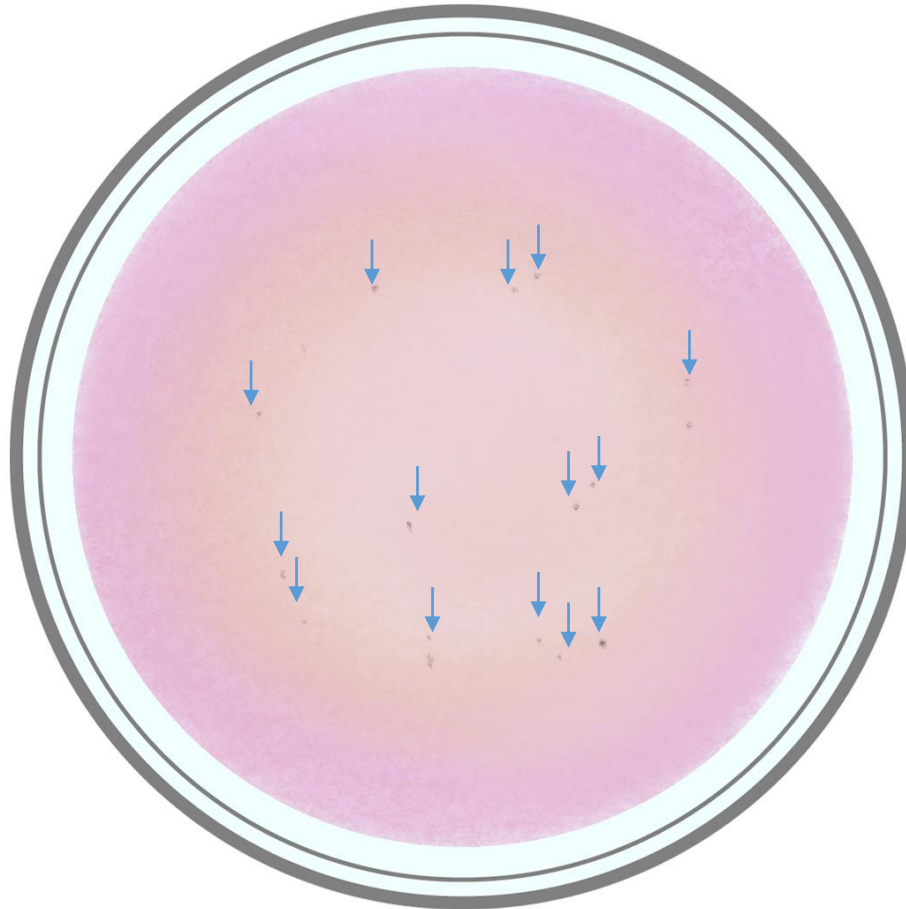


ETB


SL Interpretations



SL

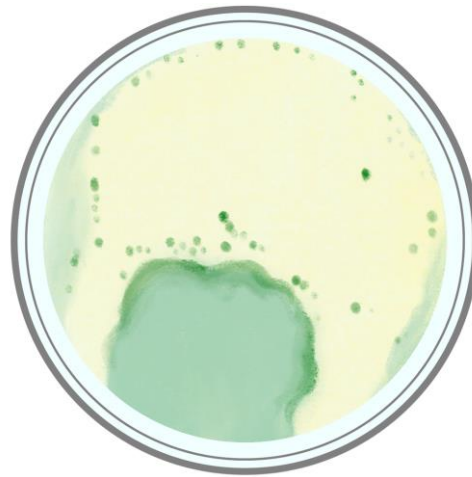


SL

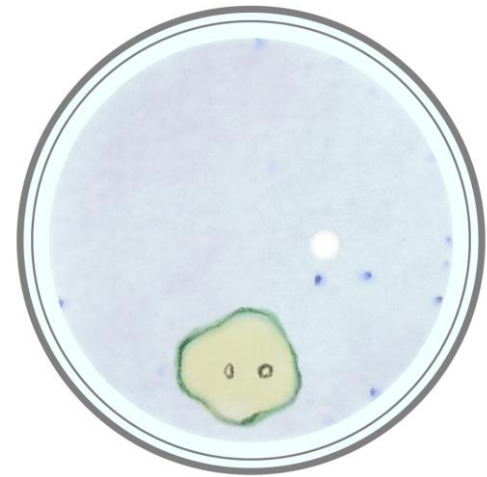


SL

SL Interpretations

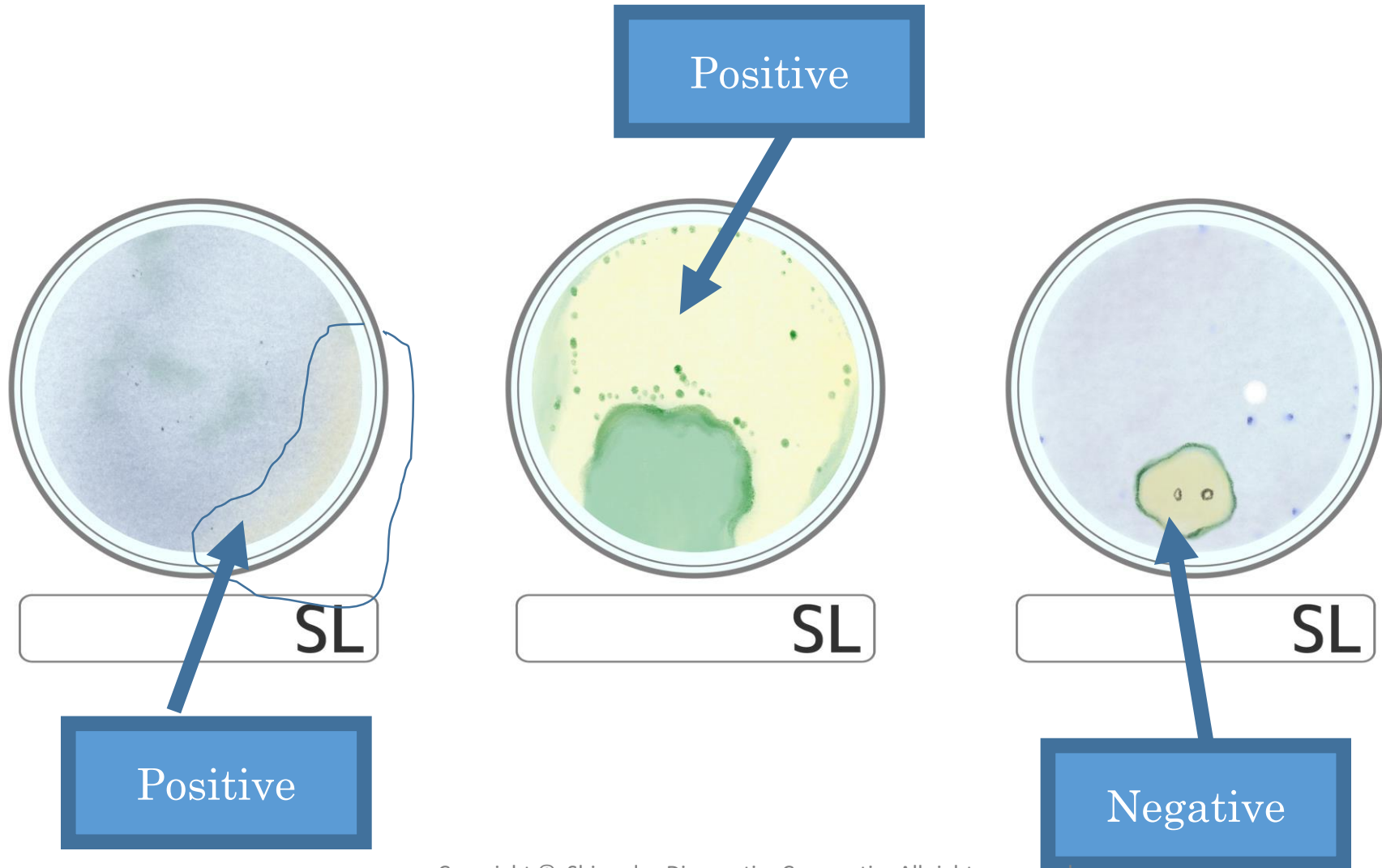


White Colonies

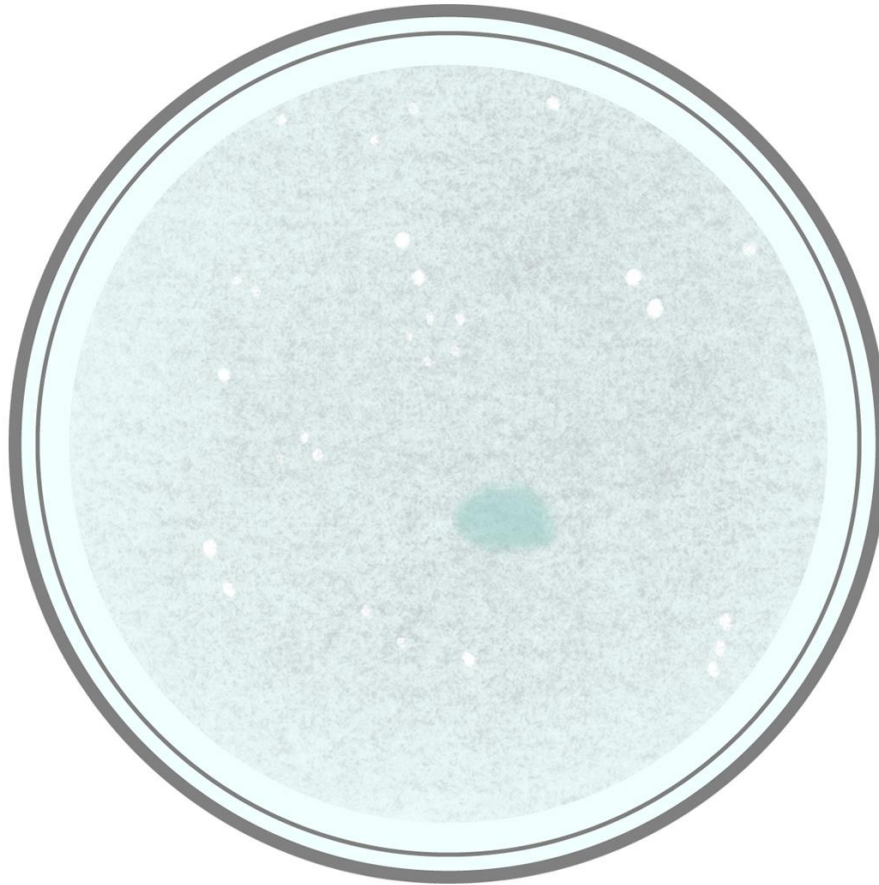

XSA


White Colonies

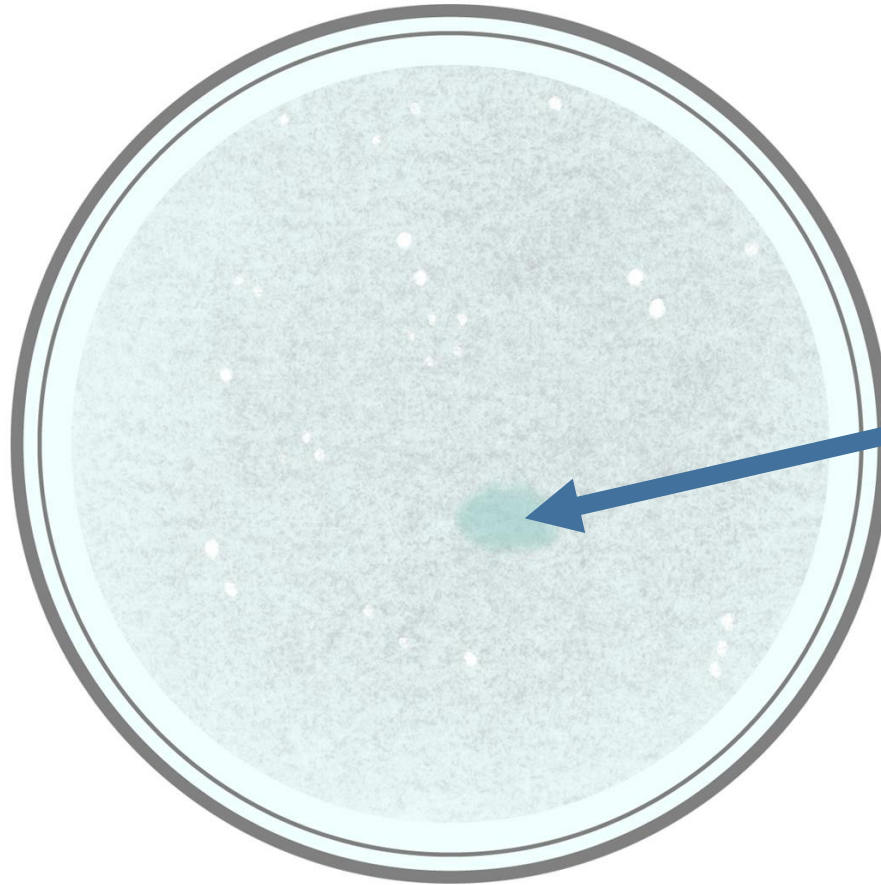

Positive


XSA

White Colonies

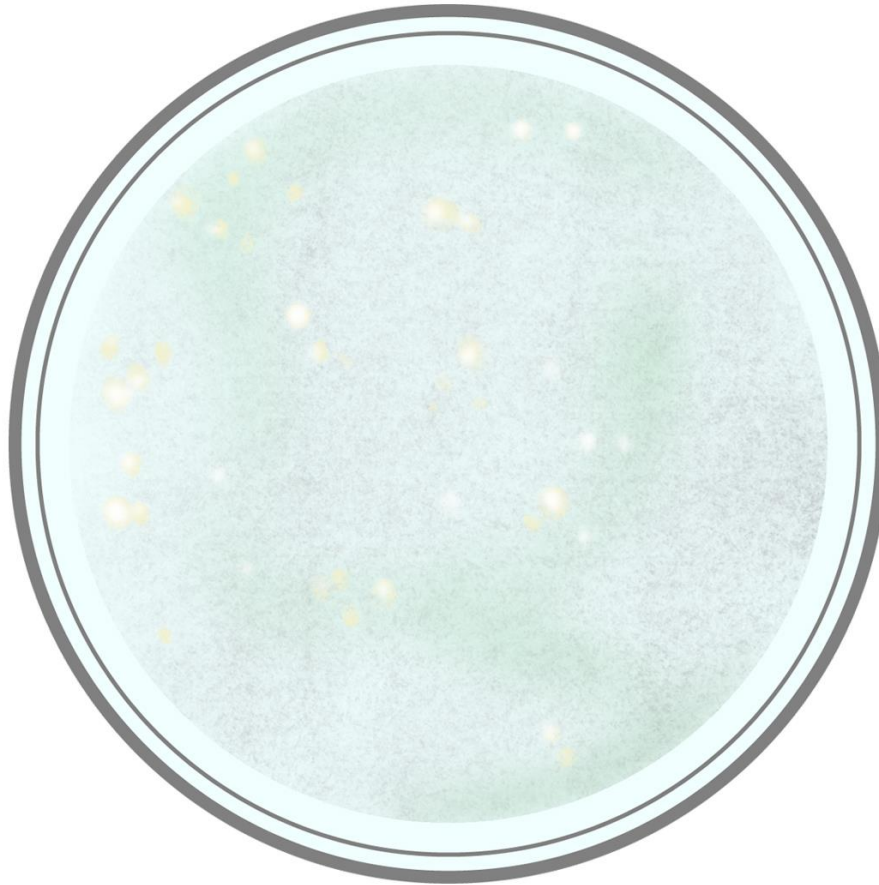

LS


White Colonies

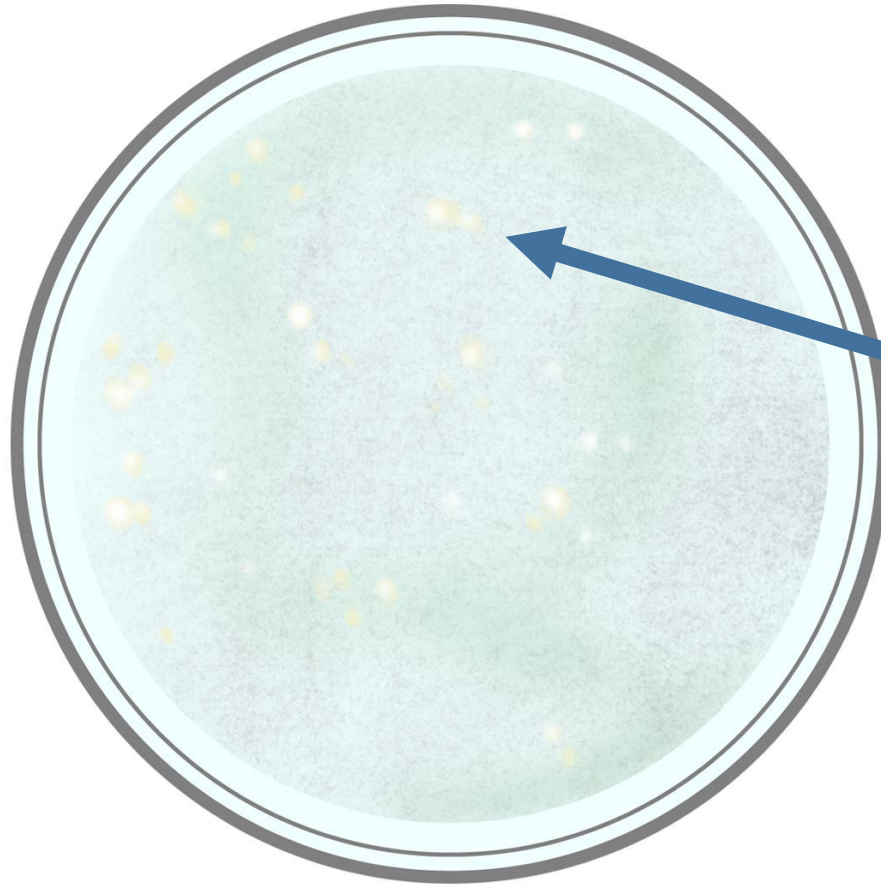

Negative


LS

White Colonies

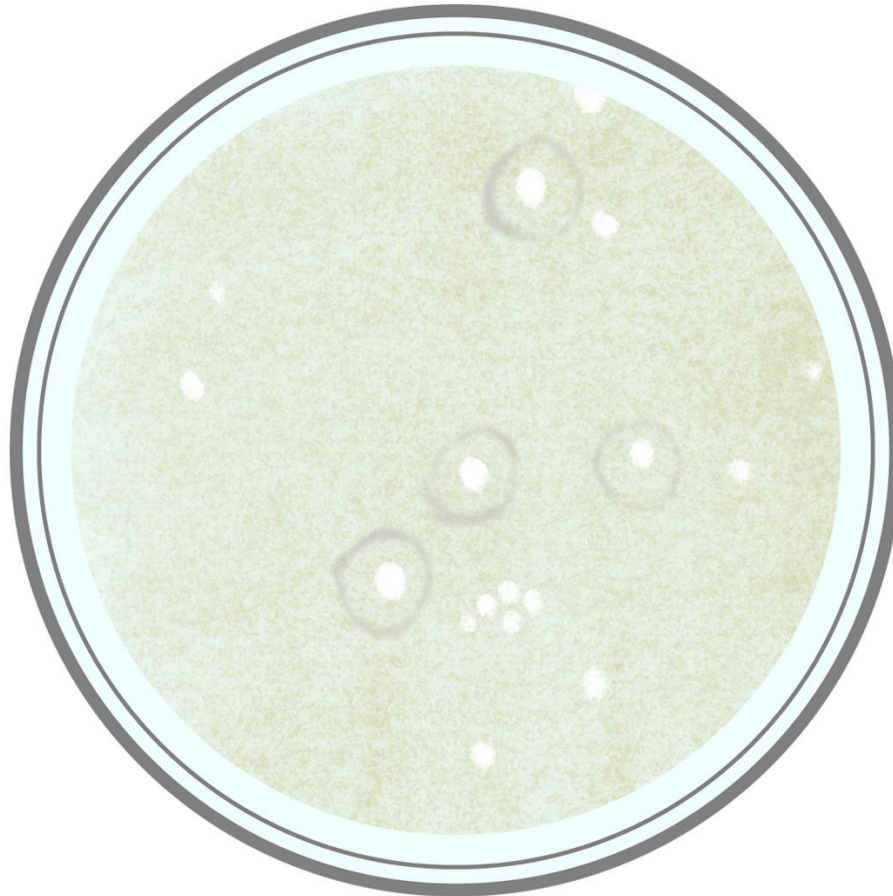

MY


White Colonies

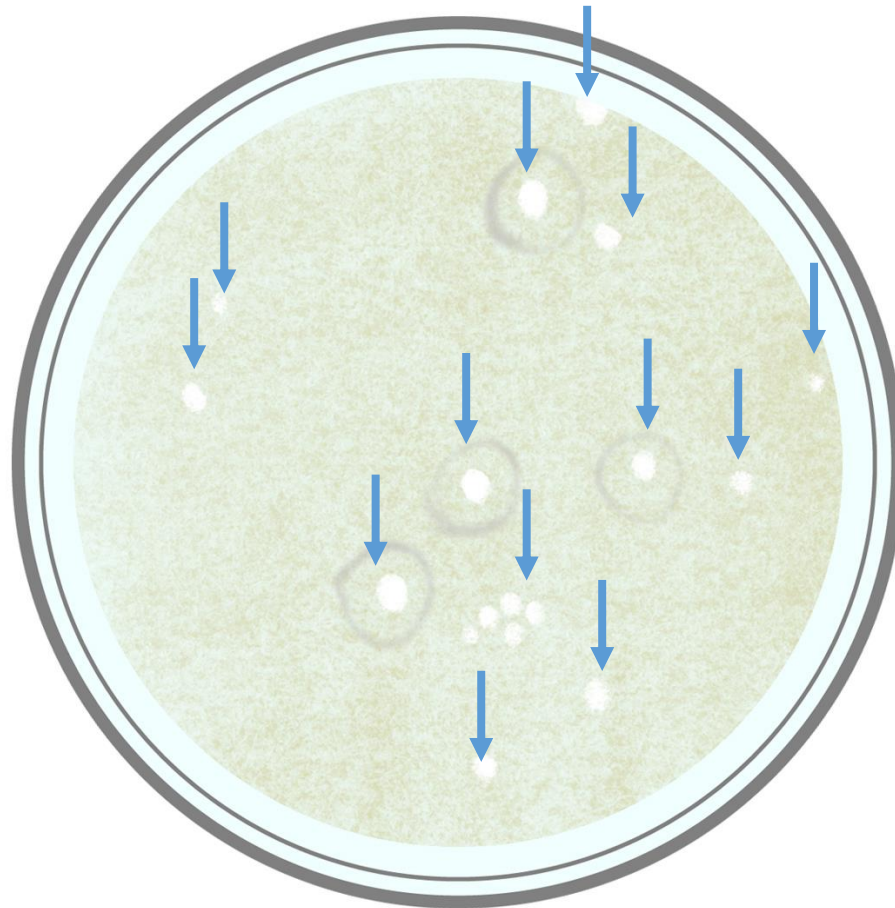

Yeasts


MY

Different Colors' Colonies

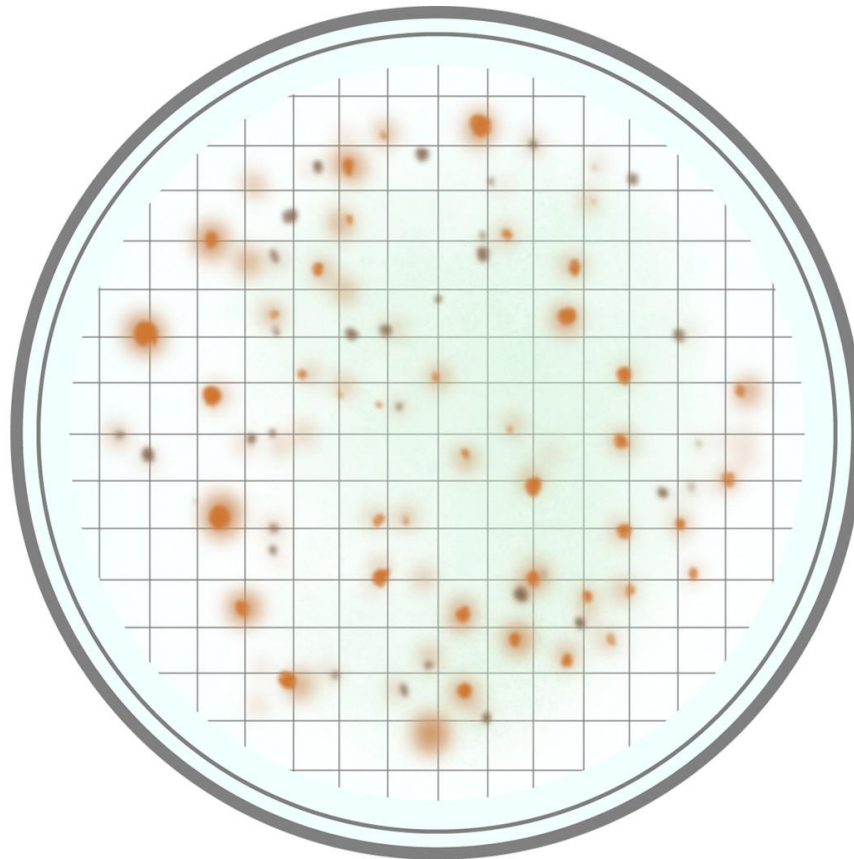

PA

Different Colors' Colonies

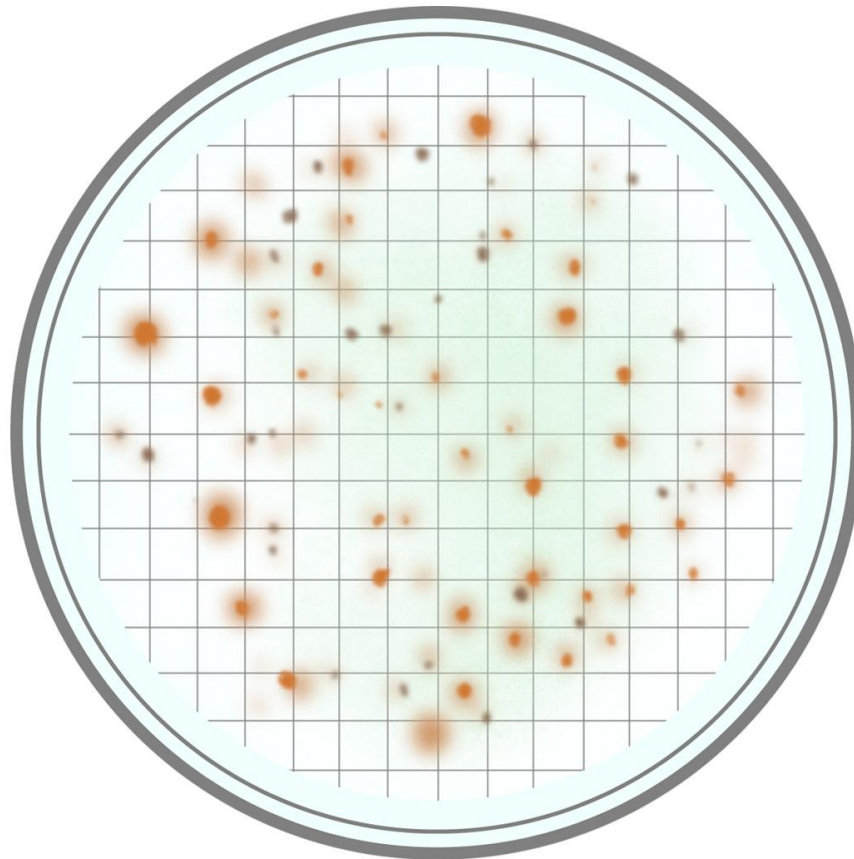

Negative


PA

Different Colors' Colonies



VP

Different Colors' Colonies

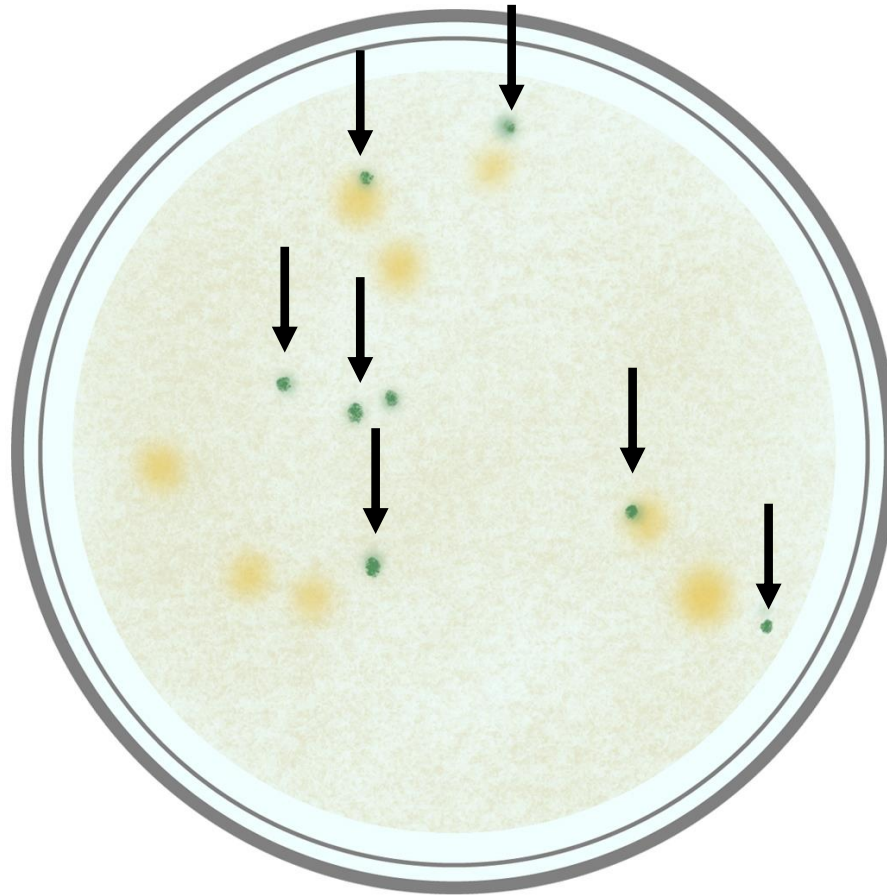


VP

Different Colors' Colonies



XSA

Different Colors' Colonies

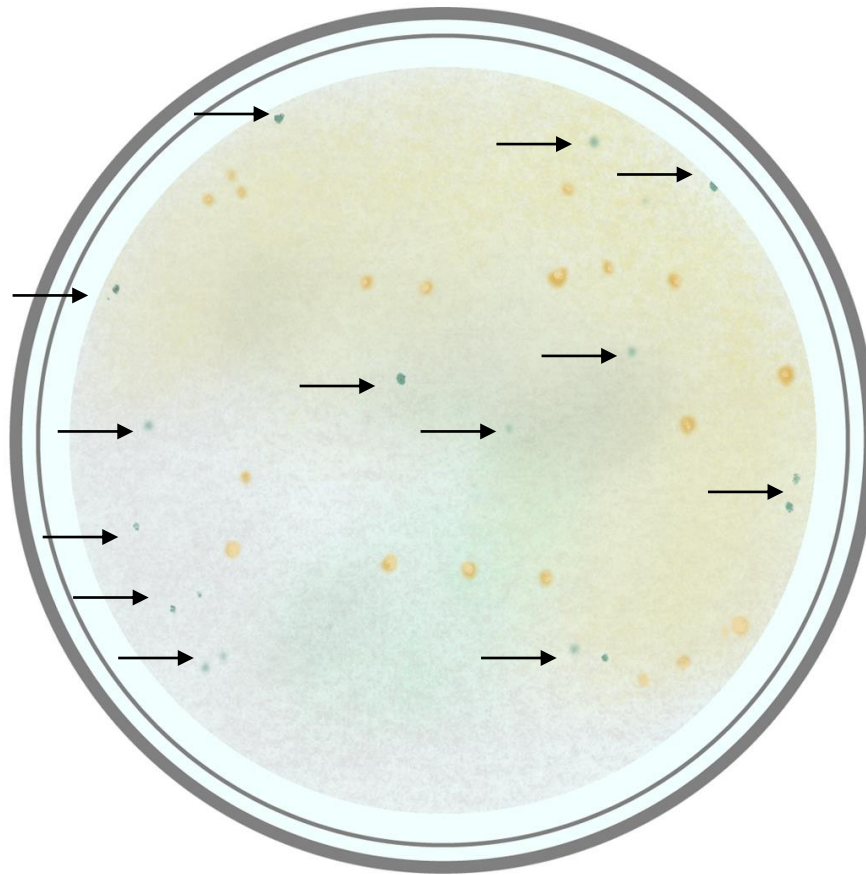

XSA