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



Compact Dry AQ Illustration Manual

Shimadzu Diagnostics Corporation




Weigh 50g solid sample

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and add 450mL Buffering Solution to the sample.





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
It is recommended to use a stomacher bag with filter to eliminate risks of carry over of tiny pieces of foodstuffs into the surface of the medium.





Homogenize this mixed sample by a blender

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Open aluminum bag, and take out a set of 4 plates.







Take off the cap of the plate

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
Write the appropriate information on the memorandum section.




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
Pipette 1ml of homogenized specimen (to be further diluted if necessary) in the middle of dry sheet of Compact Dry AQ.



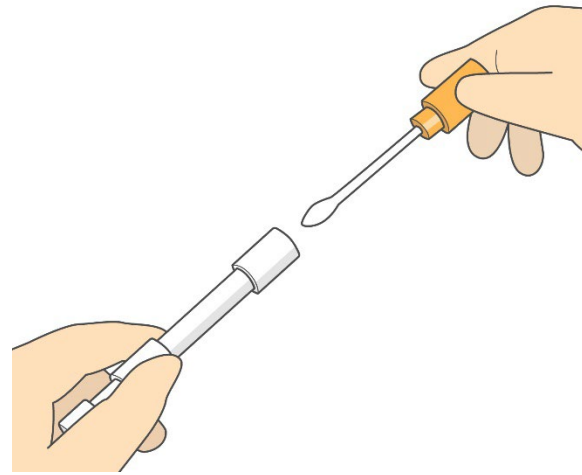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
Specimen diffuses automatically and evenly into all over the sheet (total medium of 20 cm²) to transform it into gel within seconds.




Viable count in swab
test sample

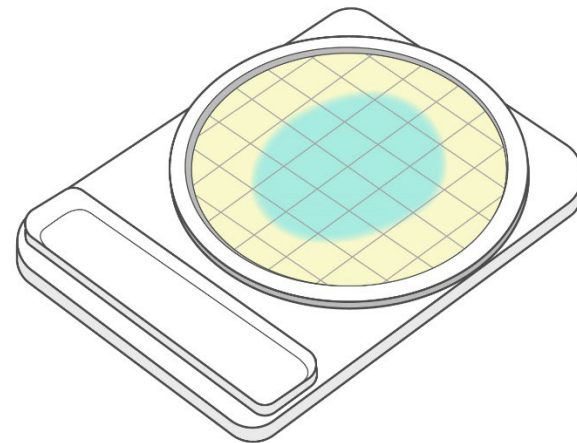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Inoculate 1 ml of
wiping solution (to be
diluted if necessary),
which is obtained from
cotton swab,



Turn over the plate capped

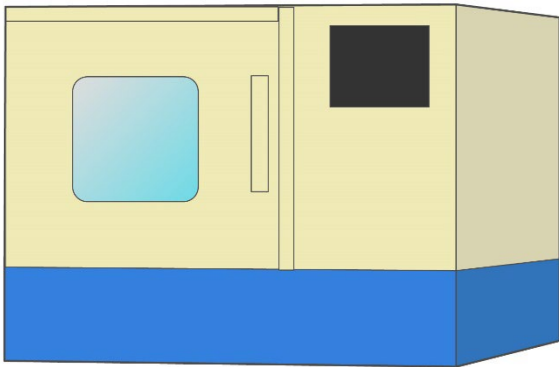

put in an incubator.

Incubate for 44 ± 4 hours at 36 ± 2 °C.

Another plate is incubated for 68 ± 4 hours at 22 ± 2 °C.